EN

PT

SV

DA

PL

NL

ValorTM 1000 Series

User Guide

1. SAFETY INFORMATION

Definition of Signal Warnings and Symbols

WARNING	For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
CAUTION	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
Attention	For important information about the product.
Note	For useful information about the product.

Warning Symbols

General hazard

Electrical shock

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

2. INSTALLATION

2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.

2.2 Connecting Power

AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the plug to a suitable electrical outlet.

Attention: Only use an AC adapter specified by OHAUS.

DC jack

Rechargeable Battery Power

Allow the battery to charge for 12 hours before using the scale on battery power. The instrument can be operated during charging, and the battery is protected against overcharging. For maximum operating time, the battery should be charged at room temperature.

2.3 Levelling the Instrument

To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.


2.4 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

3. OPERATION

3.1 Controls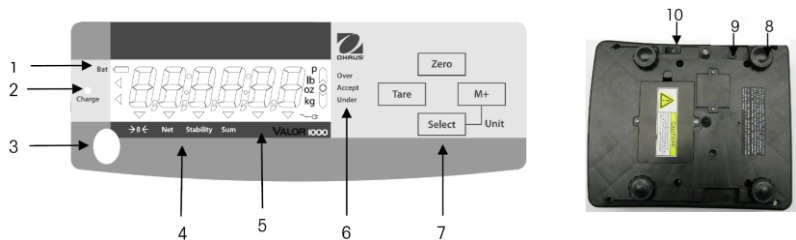

CONTROLS

Item	Description
1	Low battery indicator
2	Battery charge indicator
3	Level Bubble
4	Center of Zero, NET, Stability Sum icons
5	7-segment, 6 digit backlit LCD w/ units of measure enunciators
6	Over, Under , Accept Check weigh indicators
7	Control Buttons
8	Adjustable feet
9	Power input jack
10	On/Off Power Switch

CONTROL FUNCTIONS

Button	Action	Function
ZERO	Short Press	Sets display to zero
Enter	Menu Press	Confirms settings
TARE	Short Press	Tares weight of item on pan
	Menu Press	Decrement over and under settings
M+ Units	Short press	Stores displayed weight in accumulation memory Display accumulation data when display is at zero
	Menu Press	Increment over and under settings
	Long Press	Change unit of measure while already pressing SELECT button
SELECT Units	Short press	Shift to next digit when setting over and under values
	Long Press	Change unit of measure together with short press of M+ button

3.2 Using the Instrument

Weighing application:

1. If required, place an empty container on the pan and press Tare.
2. Add sample to the pan or container. The display shows the weight of the sample.

4. MAINTENANCE

4.1 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website www.ohaus.com to locate the Ohaus office nearest you.

5. TECHNICAL DATA

The technical data is valid under the following conditions:

Indoor use only

Operating temperature: 0°C to 40°C

Maximum Relative humidity: 80% for temperatures up to 31°C, decreasing linearity to 50% relative humidity at 40°C.

Altitude: Up to 2000m.

Power: AC adapter : 9 or 12 VDC

Pollution degree: 2

Installation category: Class III

Specifications:

Single Display Models	V11P3	V11P6	V11P15	V11P30
Dual Display Models	V11P3T/ V11P3TG*	V11P6T/ V11P6TG*	V11P15T/ V11P15TG*	V11P30T/ V11P30TG*
Capacity (kg)	3	6	15	30
Readability (g)	0.5	1	2	5
Repeatability (g)	1	2	5	10
Linearity (g)	1	2	5	10
Power requirements	9 VDC 500 mA AC adapter (supplied) Internal rechargeable sealed lead acid battery			

6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at www.ohaus.com/ce .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .

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P/N 30283480 A © 2015 Ohaus Corporation, all rights reserved.
Ohaus Corporation