**VE Series Floor Scale
Instruction Manual**

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Conformance

This product conforms to the EMC directive 89/336/EEC and the Low Voltage Directive 73/23/EEC. The complete declaration of Conformity is available from Ohaus Corporation

Disposal

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.

If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.

ISO 9001 Registration

In 1994, Ohaus Corporation, USA, was awarded a certificate of registration to ISO 9001 by Bureau Veritus Quality International (BVQI), confirming that the Ohaus quality management system is compliant with the ISO 9001 standard's requirements. On May 15, 2003, Ohaus Corporation, USA, was re-registered to the ISO 9001:2000 standard.

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INTRODUCTION

This manual contains installation, operation and maintenance instructions for the VE Series Floor Scale.

SAFETY PRECAUTIONS

Please follow these safety precautions:

- Never operate in hazardous areas.
- For VE Series models ending with S, M, or L, (Painted Steel versions) – These models may be used in dry environments only.
- For VE Series models ending with SW, MW, or LW, (Stainless Steel versions) – These models may be used in wet or dry environments.
- The foundation at the installation site must be capable of safely supporting the weight of the floor scale plus the weight of the maximum load.
- Ensure that vibrations due to machines near the installation site are kept to a minimum. This applies also when the floor scale is integrated into conveying systems.

INSTALLATION

CAUTION: THE FLOOR SCALE IS VERY HEAVY, ALWAYS USE AN APPROPRIATE HOIST TO LIFT IT FROM THE PACKAGING OR TO MOVE IT INTO THE REQUIRED POSITION.

Unpacking

1. Remove the outer packaging materials from the top and sides of the floor scale.
2. Install the two included eye bolts into opposing sides of the floor scale, see Figure 1.
3. Lift the floor scale from the packaging material, see caution notice.

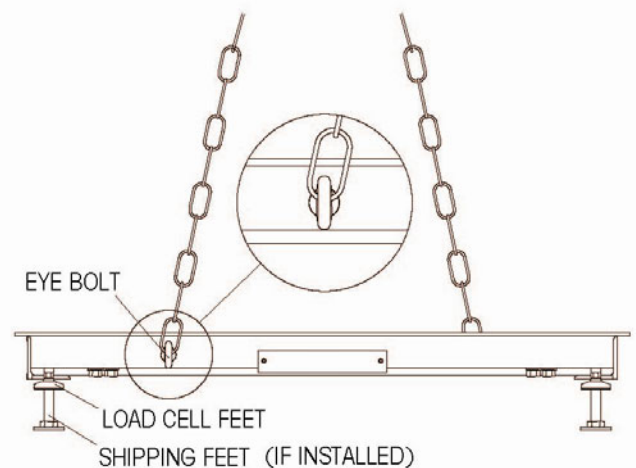
Figure 1. Hoisting Method.

Positioning

Ensure the surface is flat in the area where the scale will be positioned, especially in the region of the load cell feet. Small differences in height can be compensated by adjusting the load cell feet.

1. If an Indicator is included with the floor scale ensure that it is kept with the floor scale during positioning as there may be an attached connecting cable.
2. With the floor scale positioned over the final work area, remove the 4 shipping feet (if installed).
3. Install the 4 load cell feet (if not already installed).
4. Lower the scale on to the final work area.
5. Check that the floor scale is level; check that the 4 feet are in contact with the floor; make adjustments by rotating the nuts on the load cell feet. Use an external level.
6. Ensure that the floor scale's output cable does not become trapped or crushed during lifting or positioning.

Certifiable Installations

When used in certifiable applications at least one set of Foot Plates must be installed, (available separately). Foot Plates mount around the feet of the Floor Scale and prevent movement of the scale. Installation of Foot Plates must be done before final placement of the Floor Scale. Refer to the section on Accessories for Foot Plate installation.

Connecting an Indicator

Certain Floor Scale systems will have an Indicator already attached to the cable end. If the Indicator is not attached or if the Indicator is purchased separately follow the steps listed below to make the proper connections.

1. Route the Floor Scale cable to the position where the Indicator will be mounted.
2. Consult the installation instructions of the Indicator to properly attach the cable connections inside the Indicator,
3. The Signal names of the Floor Scale wires are shown in Table 1, Floor Scale Wiring Color Code.


CAUTION: Ensure that the cable is protected against possible damage when routed to the Indicator.

TABLE 1. FLOOR SCALE WIRING COLOR CODE

Indicator Signal Name	Floor Scale Wire Color
EXC+ (IN+)	Green
SEN+	Yellow
SIG+ (OUT+)	White
SIG- (OUT-)	Red
SEN-	Blue
EXC- (IN-)	Black
Shield (GND)	Yellow/Green

OPERATION

For setup, calibration and proper operation, refer to the instruction manual supplied with the indicator. Refer to the specifications section of the floor scale for capacity, and readability information.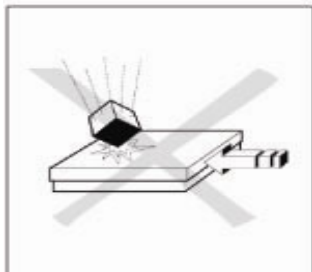

For continued best operating performance:

- Avoid falling loads, shocks and hard lateral impacts
- When weighing, all items should be centrally located on the floor scale and not overhanging the sides or the ramps.
- Verify the calibration of the floor scale system at regular intervals.

Operating Limits

Although the floor scales have extremely rugged construction, certain load limits must not be exceeded. Depending on the type of loading, the maximum safe static load, i.e. the maximum admissible loading, is shown in Table 2.

TABLE 2. MAXIMUM ADMISSIBLE LOADING.


Model	Maximum Central load ¹ (80% Surface Area)	Maximum Side load ² (30% Surface Area)	One Side Corner Load ³ (10% Surface Area)	
VE1500S/R/L	2500 kg	1500 kg	1000 kg	
VE3000R/L	3800 kg	1800 kg	1200 kg	
VE1500SW/RW/LW	2500 kg	1500 kg	1000 kg	
VE3000RW/LW	3800 kg	1800 kg	1200 kg	

MAINTENANCE

Daily Checks

- Check that the rocker feet are properly seated in the holes of each load cell.
- Check that the 4 feet are in contact with the floor.
- Check that the scale is free from debris, especially under the edges of the scale.
- Check that there is no damage to the connecting cable.


Cleaning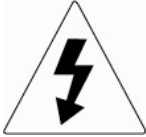

Risk of Electrical Shock


Disconnect the power source before cleaning the scale with any liquids.

Maintenance of the floor scale is limited to regular cleaning.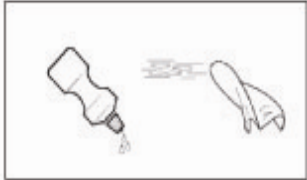

Cleaning VE-S/R/L models, (Painted Steel version) in a dry environment:

- Wipe surfaces with a damp cloth.
- Use only common household cleaning agents.
- Never use water spray or high pressure cleaners.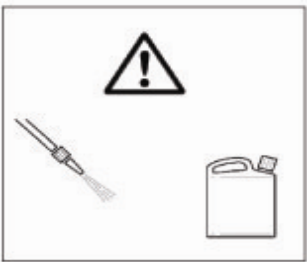

Cleaning the VE-SW/RW/LW models, (Stainless Steel version) in a wet environment:

- Water jets up to 60°C
- Use only common household cleaning agents.
- Ensure that corrosive substances are removed at regular intervals.

All models, when using optional Ramps or Foot Plates:

- Keep the air gap at the edge of the Floor Scale free of debris.

TECHNICAL DATA

Specifications

TABLE 3. FLOOR SCALE SPECIFICATIONS

Floor Scale Models	VE1500S/SW	VE1500R/RW / VE3000R/RW	VE1500L/LW / VE3000LW/LW
Dimensions	1,000 x 1,000 x 90mm	1,250 x 1,250 x 90 mm	1,500 x 1,500 x 90 mm
Capacity	1500 kg	1500 kg / 3000 kg	1500 kg / 3000 kg
Operating Temperature / Humidity	-10 °C to 40 °C / 0% to 95% Relative Humidity, Non-Condensing		
Emin	0.5 kg	0.5 kg / 1 kg	0.5 kg / 1 kg
Material, VE-S/M/L Models	Painted Carbon Steel, treaded surface, fixed top plate		
Net Weight, VE-S/M/L Models	83 kg	123 kg	170 kg
Material, VE- SW/MW/LW Models	Stainless Steel (AISI 304), sand blasted, flat surface, fixed top plate		
Net Weight, VE-SW/MW/LW Models	68 kg	101 kg	134 kg
Feet	4 x swivel design, rubber bottom, adjustable height		
Load Cells	4 x Shear Beam, Alloy Steel, OIML C3		
Load Cell Capacity Emax	1000 kg	1000 kg / 1500 kg	1000 kg / 1500 kg
Load Cell Safe Overload Limit	120% of Emax		
Load Cell Output Impedance	350 $\Omega \pm 3.5 \Omega$		
Load Cell Excitation Voltage	5V ~15V, DC/AC		
Junction Box	Rear panel access, IP65, Stainless Steel, with 5 m Cable		

DRAWINGS

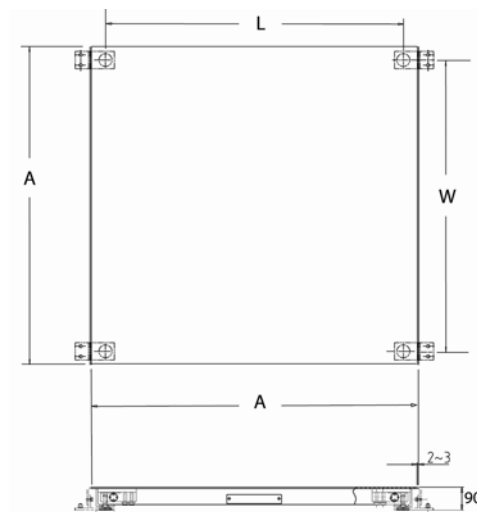
Figure 2. Floor Scale Dimensions, (Shown with optional Foot Plates)

TABLE 4. FLOOR SCALE DIMENSIONS

MODELS	DIMENSIONS (mm)		
	FLOOR SCALE PLATFORM	FOOT CENTERS	
	A	W	L
VE1500S/SW	1000	890	890
VE1500R/RW, VE3000R/RW	1250	1140	1140
VE1500L/LW, VE3000L/LW	1500	1390	1390

ACCESSORIES


TABLE 5. LIST OF ACCESSORIES

Item	Part Number
Foot Plate Kit, Painted Steel, for VE1500S/M/L	72198344
Foot Plate Kit, Stainless Steel, for VE1500SW/MW/LW	72198345
Ramp Kit, Painted Steel, 1 m, for VE1500S	72198338
Ramp Kit, Painted Steel, 1.25 m, for VE1500R, VE3000R	72198339
Ramp Kit, Painted Steel, 1.5 m, for VE1500SW	72198340
Ramp Kit, Stainless Steel, 1 m, for VE1500SW	72198341
Ramp Kit, Stainless Steel, 1.25 m, for VE1500RW, VE3000RW	72198342
Ramp Kit, Stainless Steel, 1.5 m, for VE1500LW, VE3000LW	72198343

Notes: 1. Foot Plate Kits include 2 Foot Plates, Mounting Hardware
 2. Ramp Kits include 1 Ramp, 2 Foot Plates, Mounting Hardware

VE SERIES FLOOR SCALE, INSTALLATION OF OPTIONAL FOOT PLATES AND RAMPS

Plates are required for certifiable installations. Installation of each Ramp also requires the installation of one set of Foot Plates. Foot Plates are available separately and are included with the Ramp accessory kit. Refer to Figure 3 for mounting of the Foot Plates and Ramps. Installation of Floor Plates and Ramps must be done before final placement of the Floor Scale.

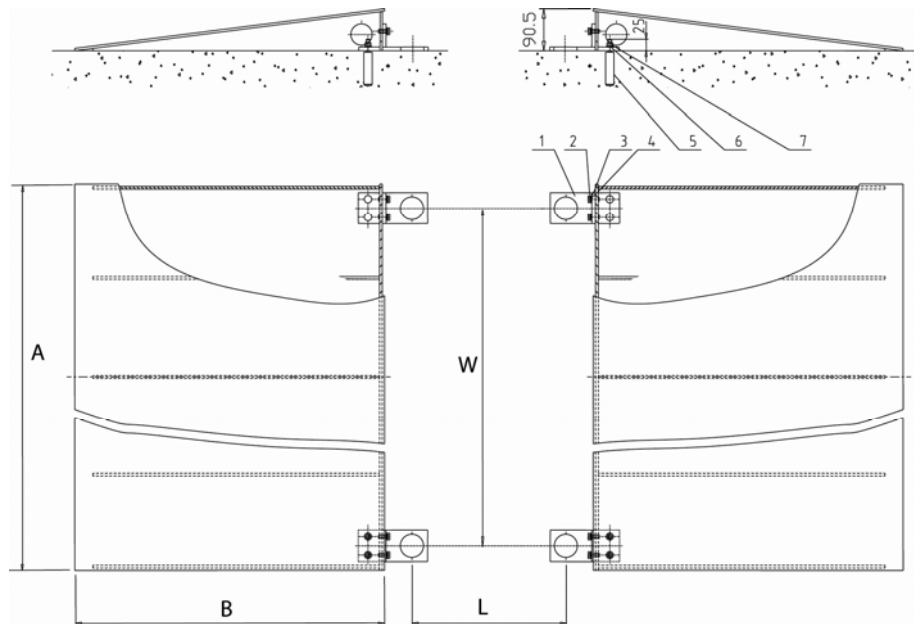LEGEND: 1. Foot Plate, 2. Bolt, 3. Spring Washer, 4. Washer, 5. Expanding Bolt, 6. washer, 7. Nut

Figure 3. Optional Foot Plate and Ramp Installation

- Notes:
1. Expanding bolts require 12.5mm diameter holes drilled 60mm deep into the floor surface.
 2. Foot Plate center to center dimensions W and L are shown in Table 4.
 3. The Floor Scale must be positioned so that the load cell feet rest within the large holes of the Foot Plates.
 4. Ensure that a free air gap exists between the sides of the Floor Scale and the edges of the Ramps.

TABLE 6. DIMENSIONS FOR FLOOR PLATE AND RAMP INSTALLATION

MODELS	DIMENSIONS (mm)			
	RAMP		FOOT PLATE CENTERS	
	A	B	W	L
Ramp for VE1500S/SW	1000	670	890	890
Ramp for VE1500R/RW & VE3000R/RW	1250	670	1140	1140
Ramp for VE1500L/LW & VE3000L/LW	1500	670	1390	1390

Warranty

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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P/N 80251484 A © Ohaus Corporation 2006, all rights reserved / todos los derechos reservados / tous droits réservés / Alle