RANGER 3000
RANGER COUNT 3000
VALOR 7000

RS232 Interface Instruction Manual

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INTRODUCTION

This RS232 interface Kit is for use with the following Ohaus products: Ranger 3000, Ranger Count 3000 and Valor 7000.

KIT CONTENTS

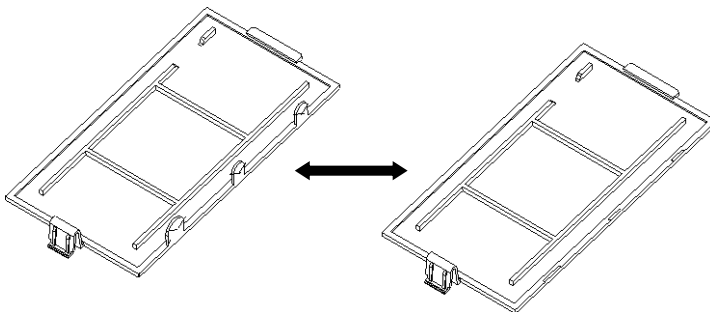
- RS232 PC Board
- Pressure Button
- Screws (4)
- Instruction Manual

INTERFACE INSTALLATION


Install the Interface module on the underside of the balance as shown.

Note: The appearance of your model may be different.

1. Remove the red interface options cover.

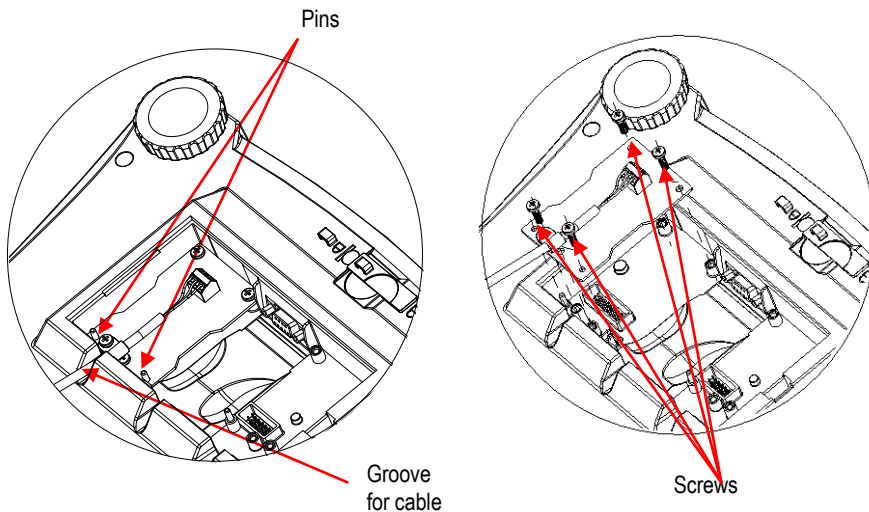2. Install the Interface PCB. Push it through the pins and secure it with the 4 screws.

SETUP

- A) The OHAUS Interface is preset to communicate using the following settings: 9600 baud, 8 bit, no parity, no handshake. Determine the RS232 parameters required for the printer or computer that is to be connected. See computer or printer documentation if assistance is required. If the parameters do not match, it will be necessary to change either the scale's settings, or the computer / printer settings.
- B) Upon installation of the interface, the scale will recognize the RS232 Interface and add relevant items to the menu. Configure the scale to the desired RS232 and printing parameters; refer to the scale User Manual for assistance in using the menus.

PRINT2

Reset
 Stable Only
 On, Off
 A.Print
 off, On.Stbl, Int, cont, Acc
 Content
 Layout
 Data.Tr
 End

COM2

Reset
 Baud Rate
 600, 1200, 2400, 4800, 9600, 19200
 Parity
 7-even, 7-odd, 7-none, 8-none
 Stop bit
 1, 2
 Handshake
 None, Xon-Xoff
 End

SETUP DEFINITIONS

PRINT / Stable - On	Balance transmits stable data only.
PRINT / Auto Print – Continuous	Balance repeatedly sends data as fast as possible.
PRINT / Auto Print – On Stable	Balance automatically outputs data upon stability.
PRINT / Auto Print – (xx) sec	Balance sends data every (xx) seconds.
PRINT / Auto Print – Off	Only sends data when PRINT is pressed.
COM / Baud, Parity, Handshake	Set parameters to match printer or computer.

CONNECTION

The RS232 Interface Kit cable terminates with a 3 pin Sub-D female connector.

Active pins: Pin 1 = Ground, Pin 2 = RXD, Pin 3 = TXD

The Interface connector will connect directly to a computer, an Ohaus printer and many other printers.

OPERATION

RS232 OUTPUT

With the Adapter installed the balance will operate in one of the following three ways:

- Auto Print = off, Stable = on or off
Pressing PRINT will send the display data.
If Stable is set to On, the balance will wait for a stable reading before sending data.
- Auto Print = not off, Stable = on or off
The balance will automatically send data based on the settings in the Menu.
If Stable is set to On, only stable values will be sent.

Data sent from the Interface is in standard ASCII format terminated with a [CRLF] (carriage return-line feed). The output format is as follows:

[weight]	9 characters (right justified)	
[space]	1 character	
[unit]	5 characters max (left justified)	
[space]	1 character	
[stability indicator]	1 character	"?" when unstable, blank when stable
[space]	1 character	
[G/N]	1 character	

Output example:

Weigh Mode Printout

	Description	Comment
11.11 kg NET	Result line	If Printx → Content → Result is ON
12.34 kg G	Gross value line	If Printx → Content → Gross is ON and a tare value is entered
11.11 kg NET	Net value line	If Printx → Content → Net is ON and a tare value is entered
1.23 kg T	Tare value line	If Printx → Content → Tare is ON and a tare value is entered
MODE: WEIGH	Mode line	If Printx → Content → Application Mode is ON
<no line printed>	Information line	If Printx → Content → Into is ON

RS232 INPUT

The balance will respond to various commands sent via the interface adapter. Terminate the following commands with [CRLF].

Commands

IP	Immediate Print of displayed weight (stable or unstable).
P	Print displayed weight (stable or unstable).
CP	Continuous Print.
SP	Print on Stability.
xS	0S: Turn off "Stable Only" menu item and allow unstable print. 1S: Turn on "Stable Only" menu item and only print stable print.
xP	Interval Print x = Print Interval (1-3600 sec), 0P turns auto print OFF.
Z	Same as pressing Zero Key.
T	Same as pressing Tare Key.
xT	Download Tare value in grams (positive values only). Sending 0T clears tare (if allowed).
PU	Print current unit: g, kg, lb, oz, lb:oz, t, c (custom).
xU	Set scale to unit x: 1=g, 2=kg, 3=lb, 4=oz, 5=lb:oz, 6=t.
xM	Set scale to mode x. M will scroll to next enabled mode.
PV	Version: print name, software revision and LFT ON (if LFT is set ON).
H x "text"	Enter Header line, where x = line number 1 to 5, "text" = header text up to 24 alphanumeric characters.
F x "text"	Enter Footer line, where x = line number 1 to 2, "text" = rear text up to 24 alphanumeric characters.
\EscR	Global reset to reset all menu settings to the original factory defaults (Escape key + 'R' Key).

AUTO-PRINT OPERATION

Once Auto-Print is activated in the menu, the balance will send data as required. To temporarily stop Auto-Printing, press the PRINT key. If there is data in the print buffer the printer will finish printing this data. A second press will resume Auto-Printing.

ACCESSORIES

For a complete listing of OHAUS printers and other accessories, contact OHAUS Corporation or visit www.OHAUS.com.

COMPLIANCE**Disposal**

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements. Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to www.ohaus.com/weee.

Thank you for your contribution to environmental protection.

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P/N 30064480B © 2013 Ohaus Corporation, all rights reserved/ tous droits réservés/ todos los derechos reservados/ Alle Rechte vorbehalten/tutti i diritti riservati