IMMB-15-02-03-22-EN-PL

Start Up Guide

Szybki start

Microbalances and Analytical Balances

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Microbalances:

Mikrowagi:

- XA 5Y.M.A
- XA 5Y.M.A.P

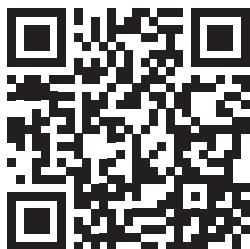
Analytical Balances:

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- XA 5Y.A

To view the complete User Manual, go to the website or scan the QR code:

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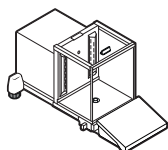
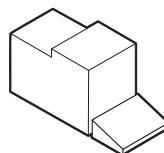
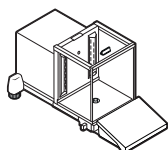
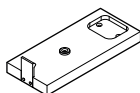
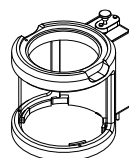
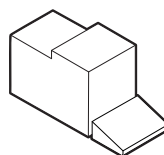
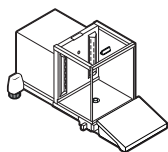
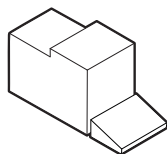
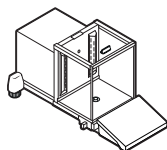
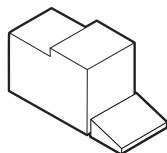
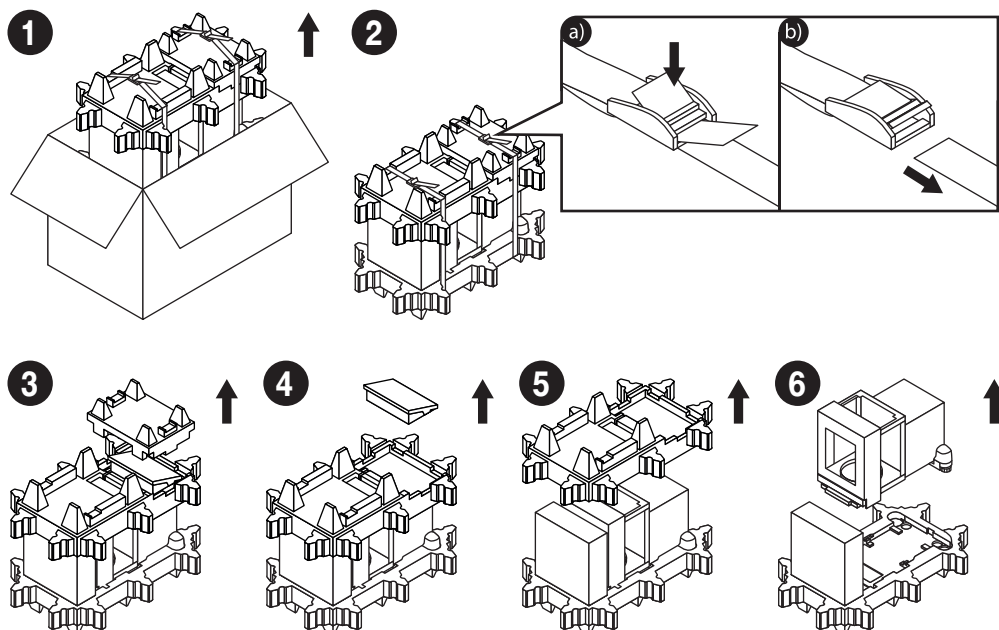
2. UNPACKING

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
Keep the packaging for warranty and service purposes.

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3. RECOMMENDATIONS AND WARNINGS

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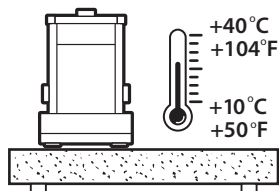
Operate the device in a room where the temperature ranges between 10 – 40 °C (50 – 104 °F) and where the relative humidity is below 80 %.

Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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
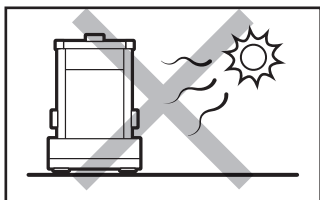
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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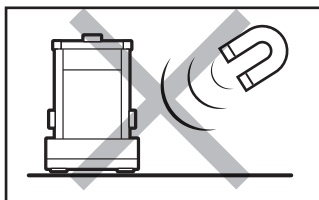
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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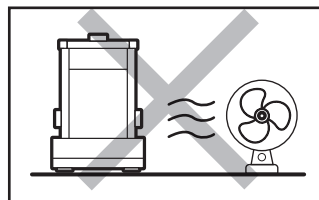
Avoid air drafts and air movements at the workstation.

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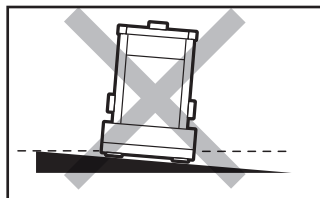
Make sure that the balance is placed on an even surface.

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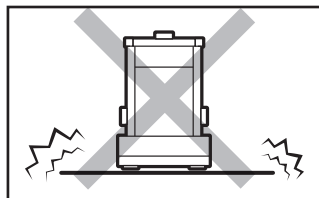Do not place the balance on an unstable ground exposed to shocks and vibrations.

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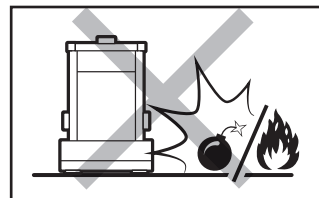Do not place the balance in a hazardous area. Do not weigh explosive materials.

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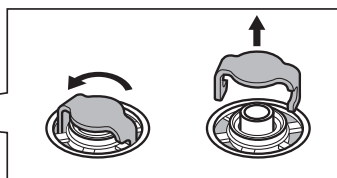
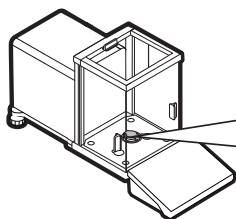
4. PRELIMINARY ACTIONS

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4.1. Transport Lock Removal

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
Remove the transport lock before assembling the balance. Keep the transport locks for warranty and service purposes.

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4.2. Grounding Spring Check

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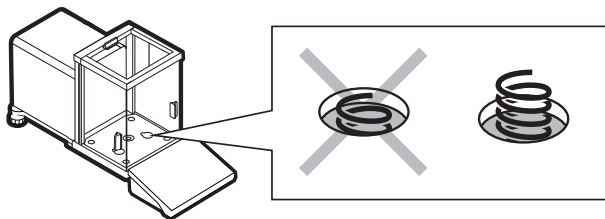Check to see if the earthing spring is placed where it is supposed to be. Make sure that the earthing spring juts slightly out of the hole.

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
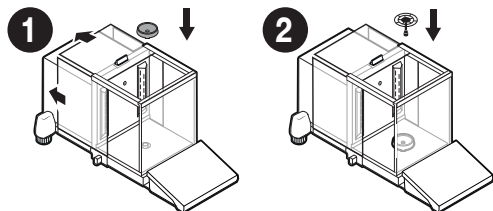
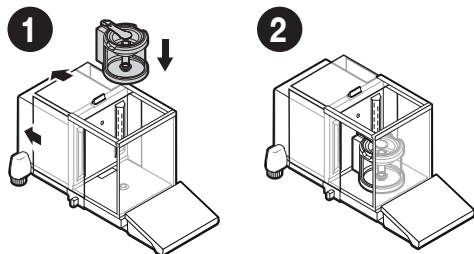
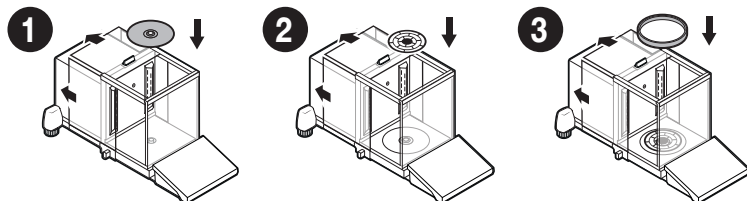
XA 5Y.A (d = 0.1 mg)

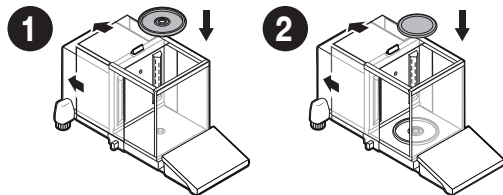6. PREPARING FOR WORK

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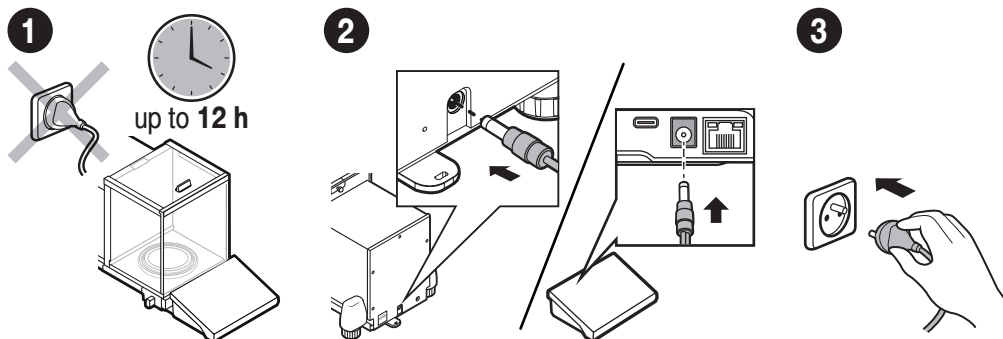
Place the device at the workstation, wait until it reaches room temperature.

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6.1. Temperature Stabilization

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Prior to switching the device on, let it reach room temperature (estimated stabilization time: up to 12 hours).

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Connect the power supply cable to the DC socket on the back of the balance or on the back of the terminal.

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Connect the power supply to the mains.

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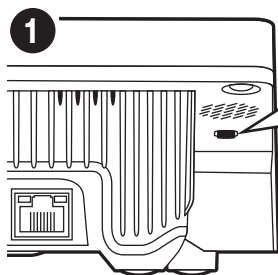
Do not connect the balance and the terminal to the power supply at the same time.

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
6.2. Switching the Balance On

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The balance switches on automatically when connected to the mains. If the balance is manually switched off using the top menu, it must be switched on using the button underneath the display housing on the left.

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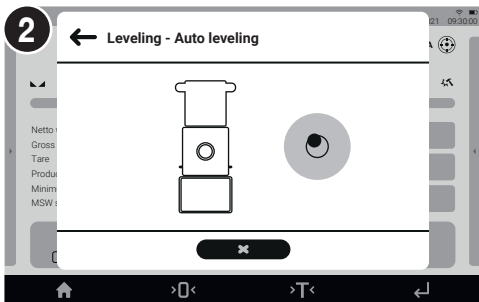6.3 Manual Levelling

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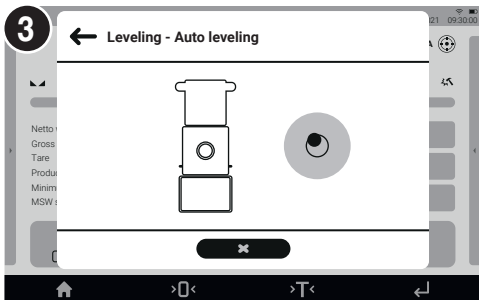
Microbalances and analytical balances are equipped with the Reflex Level System, which monitors and automatically corrects the balance level.

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

Press  to confirm the automatic levelling start. Press  to abort the process.

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
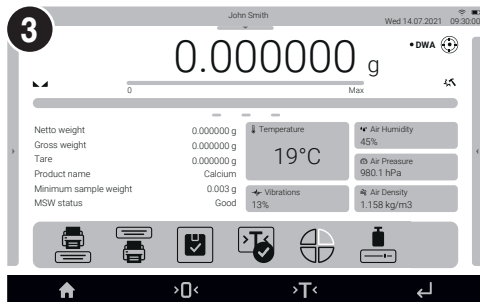
When autolevelling is complete, the home screen is displayed.

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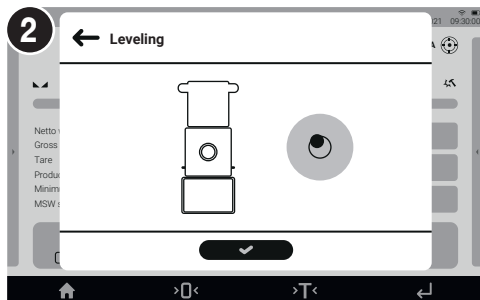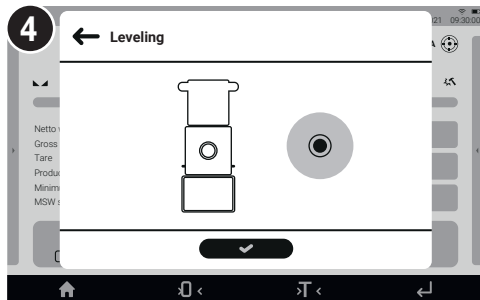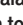The balance is levelled correctly. Press  to return to the home screen.

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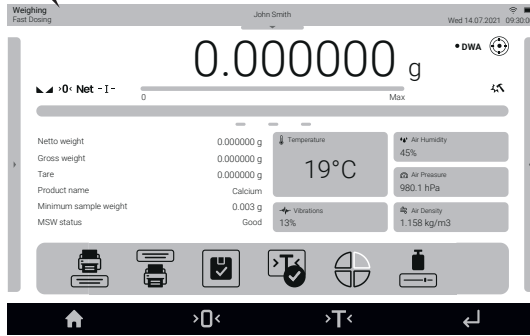
7. HOME SCREEN

EKRAN GŁÓWNY

information about the currently used working mode, logged-in operator, date, time, active connection with a computer.

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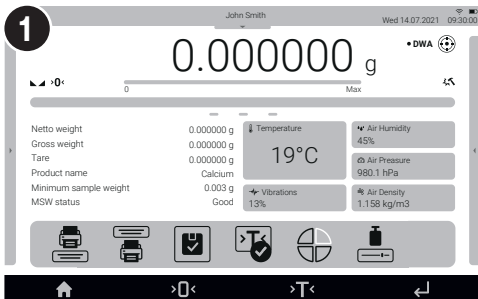
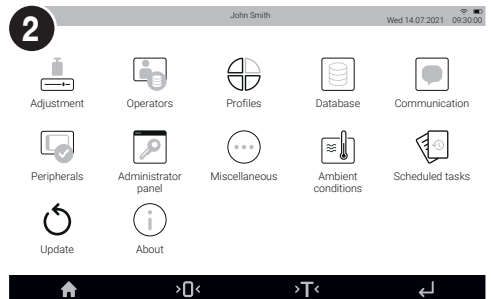
8.2. Zeroing Zerowanie

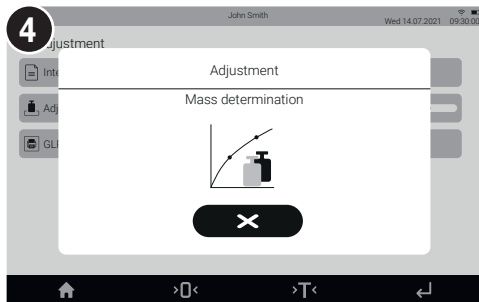
Make sure that the weighing pan is empty and press “Zeroing” button.

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
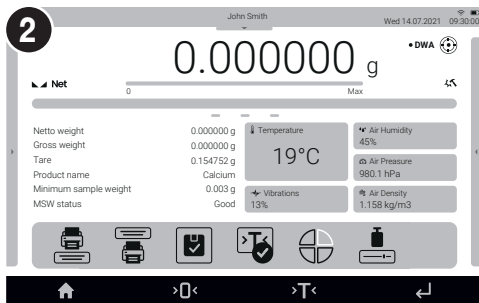
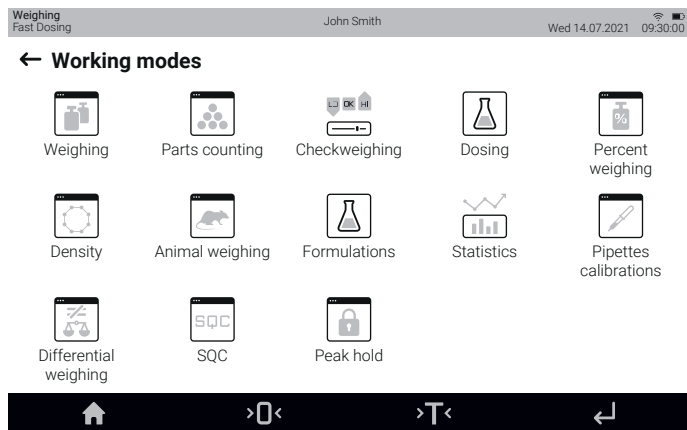
8.5. Unit Selection

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Press the current mass unit pictogram to display the list of available units.

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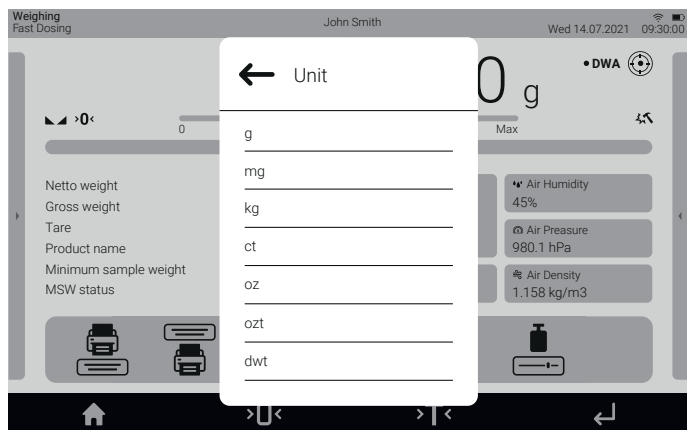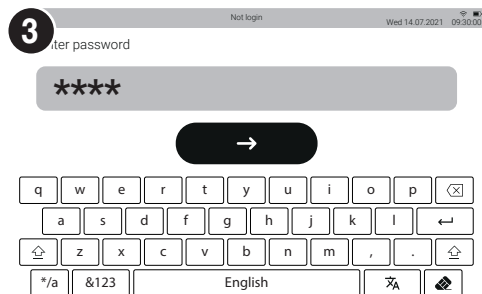
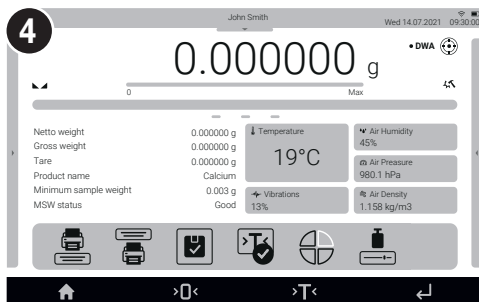
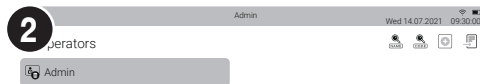
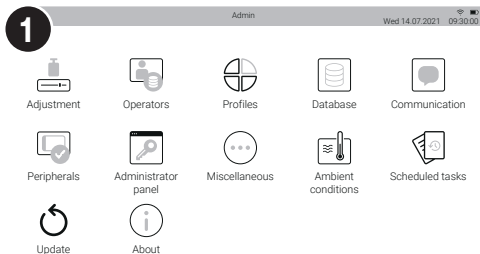
9.2. Operator Settings

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Prior to balance users setup, log in as Administrator (point 9.1).

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Hasło:

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Permissions

Assign permissions (guest, operator, advanced operator, administrator)

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Motyw

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Name and surname

Enter an operator's first and last name

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
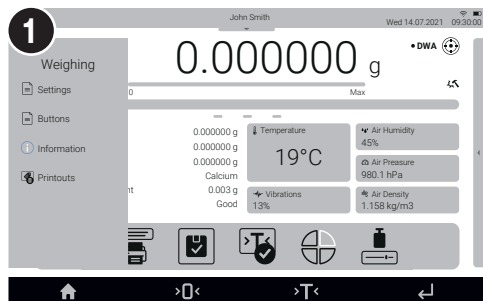
You can assign a function of your choice to each proximity sensor. In addition, you can assign various functions to the same sensor in different working modes. See the exemplary procedure for sensors configuration for the „Weighing” mode.

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
Slide out the menu.

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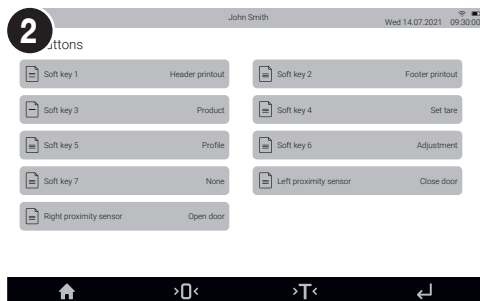Select „Buttons”.

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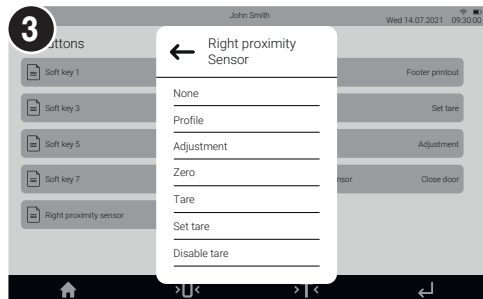
Select a respective sensor (right or left) from the list.

Select a function to be assigned to the sensor.

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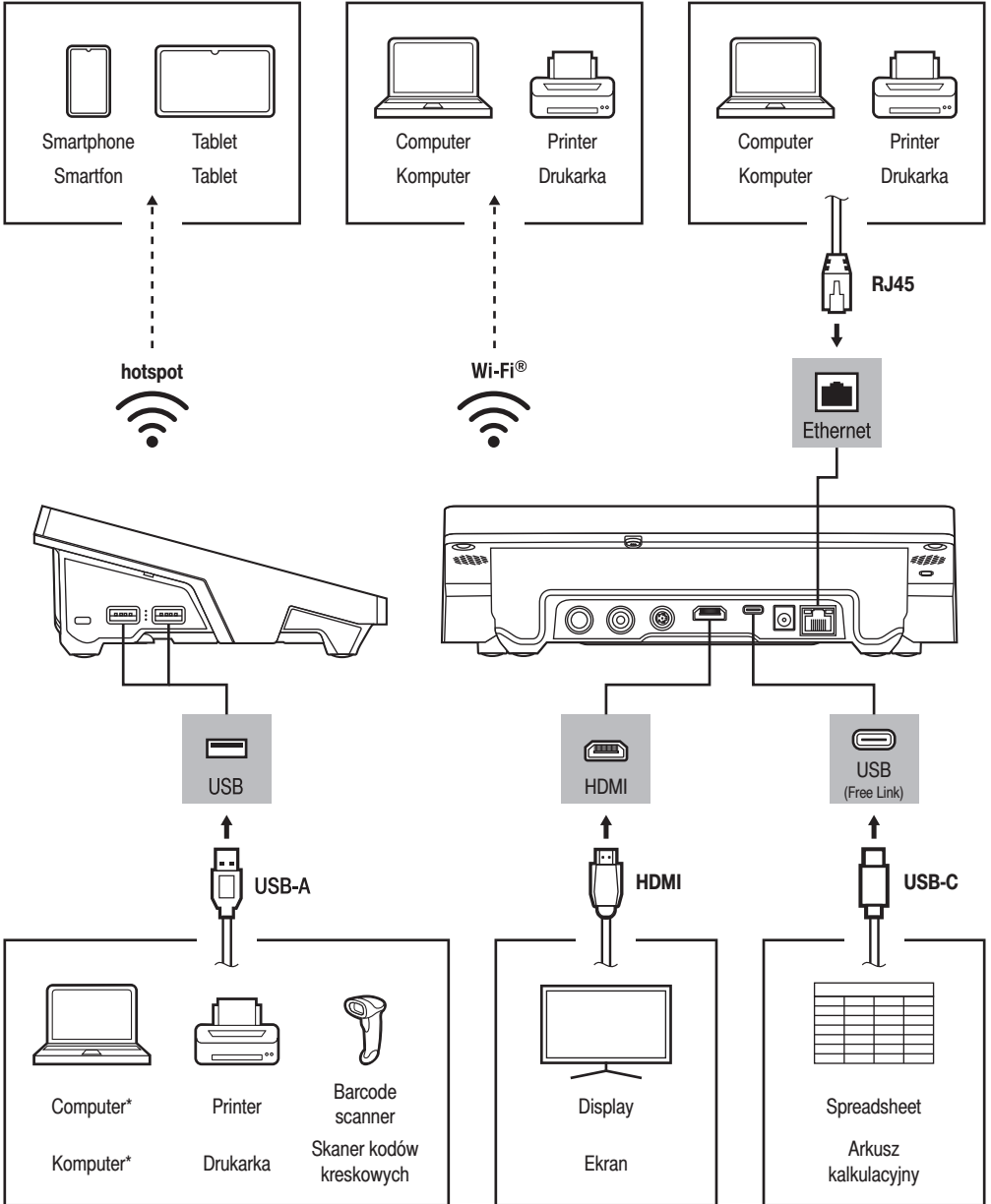Wi-Fi® is a registered trademark of Wi-Fi Alliance.

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*A USB to RS232 converter must be used for proper connection.

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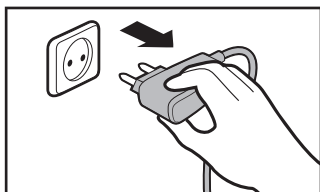
11. CLEANING AND MAINTENANCE

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
11.1. Guidelines and Precautions

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
Prior to cleaning, it is necessary to disconnect the device from the mains! Remove the plug from the outlet and disconnect all cables.

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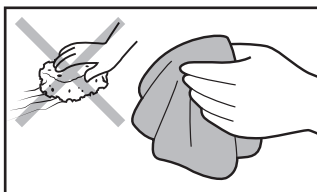Use soft cloth made of either natural or synthetic fibers exclusively. Avoid using stiff or hard fabric, as this may cause scratches on the surface!

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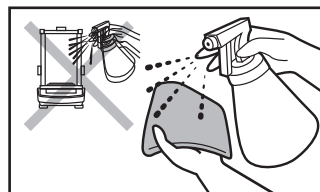Apply the cleanser to the cloth. Remember not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Use neither sharp-structured nor scrubbing-purposed substances.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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If it is necessary to disassemble the weighing chamber, follow the instructions in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either a brush or a handheld vacuum cleaner.
- Clean components can be dried using a soft cloth or a dust-free paper towel. With this, the remaining moisture will be absorbed.
- **Install all the components, providing they are completely dry.**

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
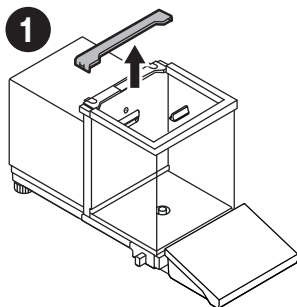
Prior to disassembling the weighing chamber elements, make sure the balance is turned off and unplugged from the mains. When disassembling the elements, use the tweezers that come standard with the instrument.

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
Grab the grips of the back pane and pull it out.

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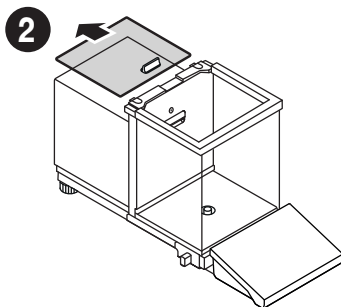Pull the top pane out.

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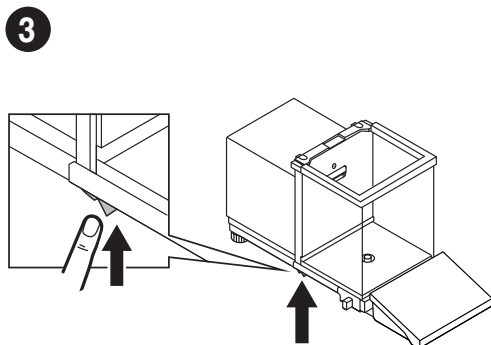Press the latch under the door to release the lock.

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