SF40A PRINTER

User Guide

EN

PT

SV

DA

PL

NL

1. SAFETY INFORMATION

Definition of Signal Warnings and Symbols

- WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
- CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
- Attention** For important information about the product.
- Note** For useful information about the product.

Warning Symbols

General hazard

Electrical shock

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for printing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

2. INSTALLATION

2.1 Connecting Power

AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

Attention: Only use an AC adapter specified by OHAUS.

2.2 Connecting the printer to the instrument

Before connecting, make sure the instrument and printer have the same RS232 parameter setting. For instrument Baud rate settings, please see the related manual.


- Attach the interconnect cable to the printer interface of the instrument and to the printer.
- Switch the instrument on.

2.3 Switching the printer on

- Connect the printer to the mains power as described in section 2.1.
- Use the main switch <I/O> to power on.
- RS232 interface only: The pilot lamp will flash until an instrument is connected.

2.4 Inserting Paper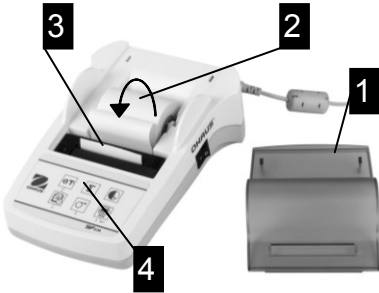

- Remove paper cover (pull on the back) (1).
- Insert the paper rotary axis through the core of the paper roll (2).
- Feed paper through the slot in the printing unit in the direction of the arrow (3).
- Press and hold until enough paper has been fed (4).
- Replace the paper cover (1).

3. OPERATION

Controls

Controls


Connections

- 1 On/Off Switch
- 2 Pilot lamp – status indicator
- 3 Control panel
- 4 Power connection
- 5 Interface connector


Control panel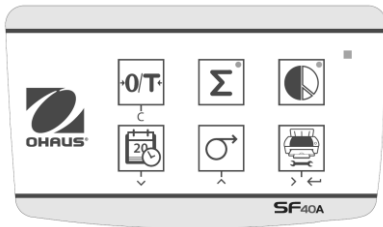

Function

Key	SF40A
	<ul style="list-style-type: none"> • Zeroing /Taring the balance • Abort – Exit
	<ul style="list-style-type: none"> • Printout of date and time • Scroll down menu option
	<ul style="list-style-type: none"> • Paper feed • Scroll up menu option
	<ul style="list-style-type: none"> • Printing of stable weight values • Opening the menu <ul style="list-style-type: none"> - select next menu option • Save settings
	<ul style="list-style-type: none"> • Totalling application
	<ul style="list-style-type: none"> • Statistics application

3.2 Menu Structure

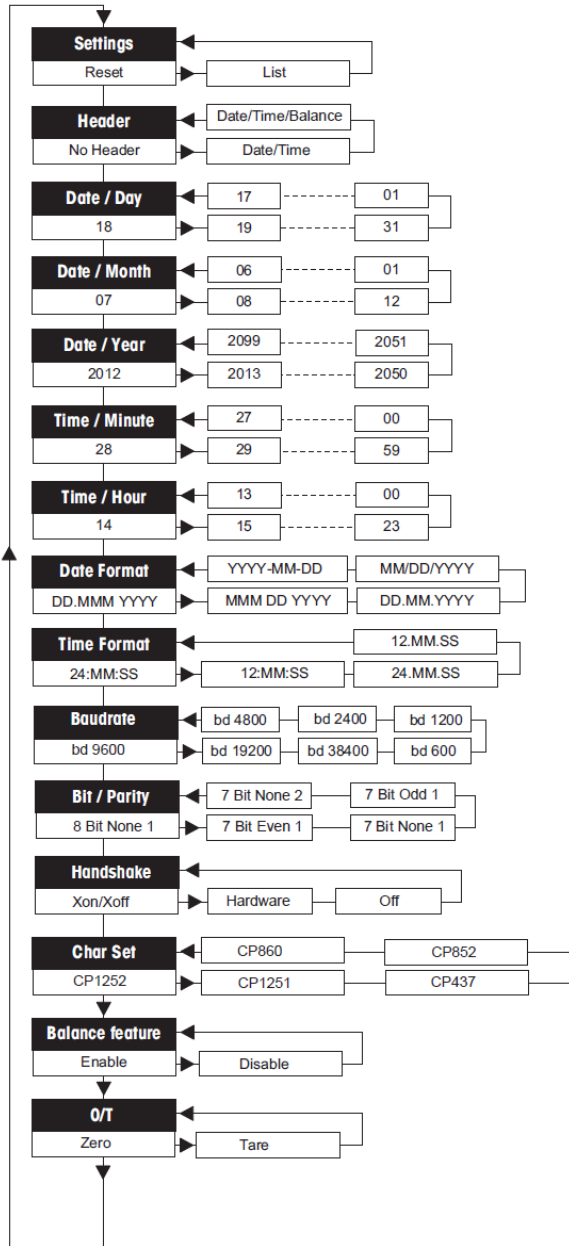3.3 Using the Instrument

Weighing application:

1. If required, place an empty container on the pan and press Tare.
2. Add sample to the pan or container. The display shows the weight of the sample.

4. MAINTENANCE

4.1 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

4.2 Troubleshooting

For technical issues contact an Authorized OHAUS Service Agent. Please visit our website www.OHAUS.com to locate the OHAUS office nearest you.

5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only.
- Operating temperature: 0°C to 40°C.
- Relative humidity: maximum relative humidity 80 % for temperatures up to 31 °C, decreasing linearly to 50 % relative humidity at 40 °C, noncondensing .
- Altitude: Up to 2000 m.
- Power input: 100-240 VAC, 1 A, 50/60 Hz", 12 VDC, 2.5 A
- Mains supply voltage fluctuations: up to ± 10% of the nominal voltage.
- Pollution degree: 2

Specifications:

Power supply	AC/DC adapter: 12 VDC, 2.5 A Primary: 100 – 240 VAC, -50 -60 Hz Voltage fluctuations: up to ±10% of the nominal voltage Secondary: 12 VDC, 2.5 A (with electronic overload protection)
Power line voltage	12 VDC
Power consumption	Max. 2.5 A

6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

	<p>This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee.</p>
--	--

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P/N 30283482 A © 2015 OHAUS Corporation, all rights reserved.
OHAUS Corporation